
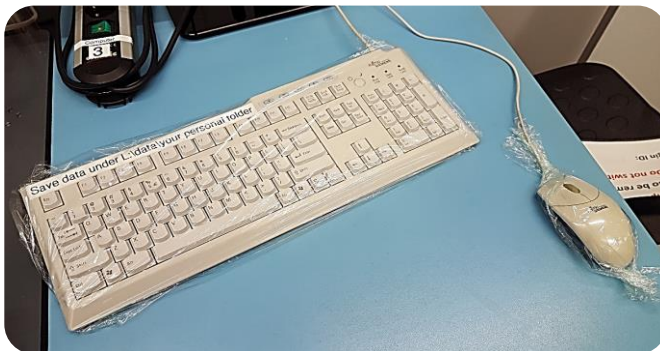


# Working at the Zeiss Imaging 3 under special safety measures

- Wear always **lab coat**, **fresh gloves**, **mask** and **safety glasses**.
- Use always **fresh plastic wrap** for this procedures!
- Cover **the control panel** before you turn on the system.
- Wrap foil around **keyboard**, **mouse**, **touchscreen** and **loosely around xyz-controller**.
- Carefully **cover** with plastic wrap foil **both oculars** and the area **above** and **below the binocular body**
- Remove **wrap** at the end of the session and **clean with 70% ETOH!**

Lab coat    fresh gloves    mask    safety glasses

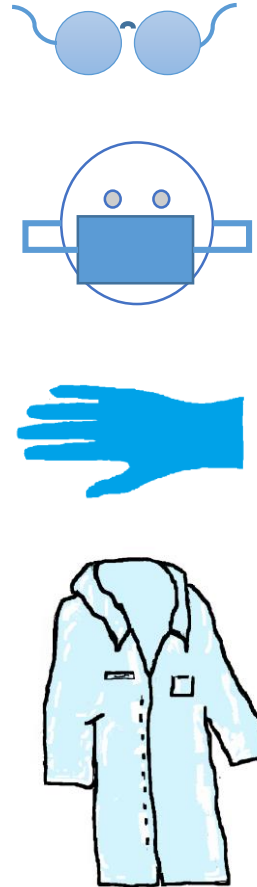

Carefully cover with plastic wrap foil **both oculars** and the **area above and below the binocular body**, use the sticky Tesafilm for fixation if needed.



**After your work** – **remove all the foils** and throw them (crumbled up) in the **yellow bin**, **clean** carefully **table surface**, **focus knobs** (at microscope body and touchscreen) and **joystick** of xyz-controller with kitchen tissue wetted **with 70% ETOH** (still wear your gloves). **Take off gloves and lab coat only when leaving LIC rooms at the main exit!**



*Please keep to these guidelines to ensure a safe working environment for all - thank you very much!*

your LIC team