
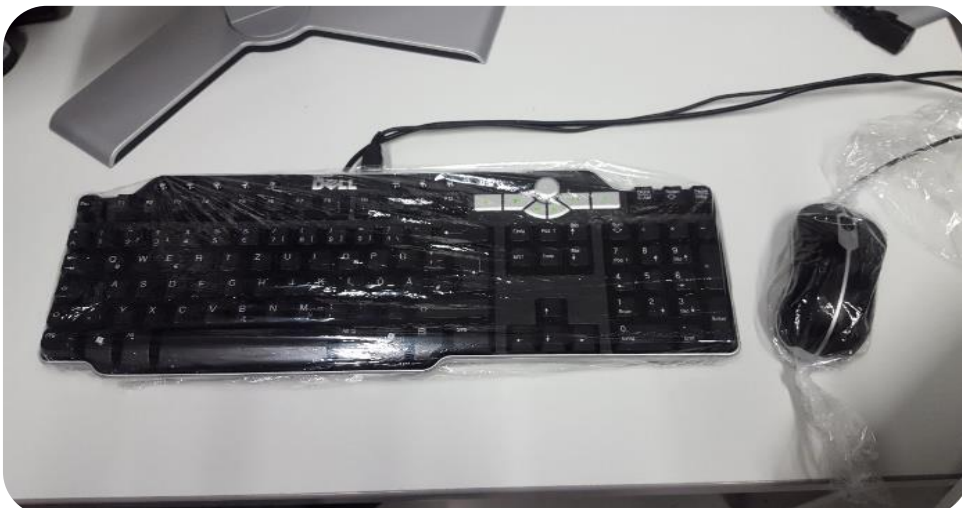
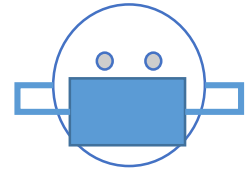



Working at the Zeiss Lightsheet under special safety measures

- Wear always **lab coat, fresh gloves, mask** and **safety glasses**.
- Use always **fresh plastic wrap** for this procedures!
- Wrap foil around **keyboard** and **mouse**.
- Remove **wrap at the end of the session** and **clean with 70% ETOH!**
- Take off gloves and lab coat only when leaving LIC rooms at the main exit!

Lab coat
fresh gloves
mask
safety glasses

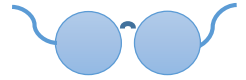

After your work – **remove all the foils** and throw them (crumbled up) in the **yellow bin**, **clean** carefully **table surface**, **focus knobs** and **joystick** of xyz-controller with kitchen tissue wetted with **70% ETOH** (still wear your gloves).
Take off gloves and lab coat only when leaving LIC rooms at the main exit!

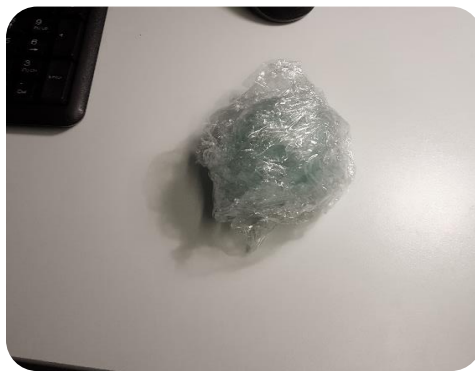

Please keep to these guidelines to ensure a safe working environment for all - thank you very much! your LIC team